

# Allergeneye™ Quick: Rapid 10 minute Surface and Rinse Water Test for Allergens

## Invisible Allergens, Visible Results

Food allergen contamination poses serious risks to consumer safety and your business reputation. Even microscopic residues left on manufacturing equipment can trigger life-threatening reactions. Traditional cleaning methods alone aren't enough – you need proof that your equipment is truly allergen-free.

## Allergeneye™ Quick: Swab or Rinse. 10 Minutes. Complete Confidence.

The Allergeneye Quick from Kikkoman is a rapid test kit designed specifically for post-cleaning allergen verification in food manufacturing environments. Whether testing equipment surfaces or rinse water after COP/CIP procedures, this innovative solution delivers reliable results in just 10 minutes.



### Key Features:

- **Results available 8 major allergens**
- **Dual Testing Capability**  
Equipment surface and rinse water analysis in one versatile kit.
- **High Sensitivity**  
Detects allergen proteins at 0.2ppm
- **Rapid Results**  
Clear determination in 10 minutes
- **Room Temperature Storage**  
Convenient storage and handling

### Allergeneye™ Quick tests are available for:

- Egg
- Milk (Casein)
- Wheat
- Buckwheat
- Peanut
- Crustacean
- Soybean
- Sesame

Learn more about  
Allergeneye™ Quick



## Sample Results in 10 minutes


### Equipment Swab Test



### Rinse Water Test



1

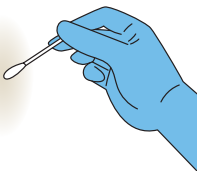


Swab the test area (10cm x 10cm) with moistened swab.




Collect rinse water after cleaning equipment and tools that are subjected to testing.

2

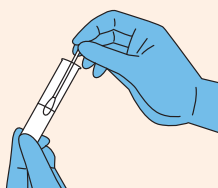


Put the swab in 10mL buffer solution and stir it thoroughly.

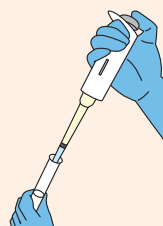


Add phosphate buffered saline to the rinse water to make a 10-fold diluted test specimen.

3

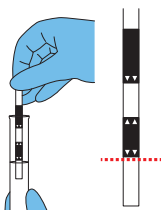

Immerse the strip up to DIP line for 3 seconds and then remove it.

4

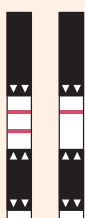

After 10 minutes, results are indicated.  
Left : positive  
Right: negative

## Allergeneye™ Quick

Measurement Time: 10 minutes

Detection Limit = 0.2ppm

Storage Condition : 1-30°C (Room Temperature)

Kit Name	Product Code	Test Target	Label Color	Tests
Allergeneye™ Quick for Egg	61962	Ovalbumin	Yellow	20
Allergeneye™ Quick for Milk (Casein)	61963	αs1-casein	Blue	
Allergeneye™ Quick for Wheat	61964	Gliadin	Red	
Allergeneye™ Quick for Buckwheat	61965	24 kDa protein	Green	
Allergeneye™ Quick for Peanut	61966	Ara h 1	Brown	
Allergeneye™ Quick for Crustacean	61967	Tropomyosin	Orange	
Allergeneye™ Quick for Soy	61968	7S globulin	Light blue	
Allergeneye™ Quick for Sesame	61969	11S globulin	Black	

\* Wiping tools (swabs, buffer solution, container) are not included in this kit.

Commercially-available swabbing equipment and dilution-purpose phosphate buffered saline for microorganism tests can be used with this kit. However, such equipment and phosphate buffered saline must not contain proteins such as peptone medium or allergens.

**kikkoman**

Kikkoman Biochemifa Company

2-1-1 Nishi-Shinbashi, Minato-ku, Tokyo 105-0003 JAPAN

info\_biochemifa@mail.kikkoman.co.jp

<https://biochemifa.kikkoman.com/e/kit/>

\*The information contained in this catalog may change without notice.

©2025 Kikkoman Corp. (5112Y251101)